BACTERANIOS SF IP sterile concentrate

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- **Biocidal product with broad spectrum**
- **Synergistic combination of 2 antimicrobial active ingredients**
- **Formaldehyde free**
- **Activity in compliance with European standards**
- **Double-bagged packaged**
- **Gamma irradiated**

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INDICATIONS

Cleaning and disinfection of floors, walls, surfaces and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

Glutaraldehyde (CAS N°. 111-30-8 : 50 mg/g), didecyldimethylammonium chloride (CAS N°. 7173-51-5 : 120 mg/g).

MICROBIOLOGICAL PROPERTIES

Bactericide in 5 min, 0.5%, 20°C (EN 1276, EN 13697).
Yeasticide in 15 min, 0.5%, 20°C (EN 1650, EN 13697).
Fungicide in 30 min, 2%, 20°C (EN 1650, EN 13697).
Active in 5 min 0.5% 20°C against BVDV (surrogate of HCV) and PRV (surrogate of HBV) .

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...): see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Liquid colourless solution
- Soluble in all proportions in water
- Density at 20°C : approx. 1
- pH of product at 20°C : approx. 4
- Non-corrosive towards metals (non oxidising).

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INSTRUCTIONS FOR USE

Concentrated solution.

Remove the first package before entering into the clean room. Remove the second package on first use.

Use at dilution from 0,5% (5ml/l) to 2% (20ml/l), i.e. 1 to 4 bottles for 12l of cold water, according to the required antimicrobial activity, respecting the indicated contact time. Apply a sufficient amount (+/- 30 ml/m²).

If necessary, remove the excess after the contact time. For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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PRECAUTIONS FOR USE

Dangerous. Follow the instructions for use (in accordance with Directive 99/45/EC and its adapted versions).

Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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