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- *Biocidal product with broad spectrum*
- *Synergistic combination of 2 antimicrobial active ingredients*
- *Aldehyde free*
- *Activity in compliance with European standards*
- *Double-bagged packaged*
- *Gamma irradiated*

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INDICATIONS

Cleaning and disinfection of floors, surfaces, walls and equipment in the Pharmaceutical Industry for clean rooms (A and B grades).

COMPOSITION

N-[3-aminopropyl]-N-dodecylpropan-1,3-diamine - N° CAS 2372-82-9 (51 mg/g), Didecyldimethylammonium chloride (N° CAS 7173-51-5 (25 mg/g).

MICROBIOLOGICAL PROPERTIES

*Bactericide in 5 min, 0,5% 20°C (EN 1276, EN 13697).
Yeasticide in 5 min, 0,25%, 20°C (EN 1650).
Active in 5 min, 0,25%, 20°C against HIV-1, PRV (surrogate of HBV), BVDV (surrogate of HCV) and H1N1.*

Biocidal activity on field surfaces and strains (bacteria, yeasts, moulds) according to EN 13697.

Other data concerning the product (material compatibilities...) : see scientific file.

For each batch, Certificate of gamma irradiation, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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PHYSICO-CHEMICAL DATA

- Density at +20°C: approx. 1
- pH of pure product at 20°C: > 11
- Non corrosive (non oxidising)

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INSTRUCTIONS FOR USE

Concentrated solution.

Remove the first package before entering into the clean room. Remove the second package on first use. Use at dilution of 0,5% (5ml/l), i.e. 1 flask for 12L of cold water, respecting the indicated contact time, for the required antimicrobial activity. Apply in sufficient amount (+/- 30 ml/m²). If necessary remove the surplus after the contact time. For frequency of use and cleaning of the application equipment, please refer to the hygiene plan in place.

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PRECAUTIONS FOR USE

Dangerous - follow the instructions for use (in accordance with Directive 99/45/EC and its adapted versions). Use the biocides with precaution. Before use, read the label and the product information.

Store between +5°C and +35°C.

Biocidal product designed for surface disinfection (Group 1 - PT2) - Reserved for professional use.

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